

Food Allergen Test Kits

Reveal® 3-D

Gliadin



Intended Use

Reveal 3-D Gliadin is intended for the qualitative analysis of gliadin residue in clean-in-place rinses, and on environmental surfaces.

The Test

Reveal 3-D Gliadin is a unique hand-held device with a 3-line readout: a control line confirms the method has been performed successfully and two other lines differentiate nondetectable, low and high levels of gliadin.

The Procedure

Rinse and environmental samples:

1. Add 0.25 mL of sample or environmental swab to an extraction buffer bottle.
2. Cap and shake for 1 minute.
3. Remove cap and fill with liquid from bottle.
4. Dip the Reveal 3-D device into the liquid in the lid.
5. Leave the cavity saturated until liquid is seen running into the test window.
6. Place device in a flat surface and allow test to develop for 5 minutes.

See package insert for complete instructions.

Product Specifications

Sensitivity:	Screens samples at 5 ppm gliadin (10 ppm gluten)
Testing time:	5 minutes
Storage:	Refrigerated 2–8°C
Tests per kit:	20

Results



Ordering Information

Prod.#	Product description
8519	Reveal 3-D Gliadin