

# Reveal 3-D

## For Total Milk Allergen



### Intended Use

Reveal 3-D for Total Milk Allergen is intended for the qualitative analysis of milk residue (casein or whey) in liquid products (e.g., juices and sorbet), clean-in-place rinses, and on environmental surfaces.

### The Test

Reveal 3-D for Total Milk Allergen is a unique hand-held device with a 3-line readout: a control line confirms the method has been performed successfully and two other lines differentiate nondetectable, low and high levels of milk.

### The Procedure

Rinse and liquid products:

1. Add Type 8 buffer sachet to sample tube.
2. Add 1 mL of sample or environmental swab to a sample tube.
3. Cap and shake for 1 minute.
4. Remove cap and fill with liquid from tube.
5. Dip the Reveal 3-D device into the liquid in the lid.
6. Leave the cavity saturated until liquid is seen running into the test window.
7. Place device in a flat surface and allow test to develop for 5 minutes.

See package insert for complete instructions.

### Product Specifications

Sensitivity:	Screens samples at 5 ppm total milk
Testing time:	5 minutes
Storage:	Refrigerated (2–8°C)
Tests per kit:	10

### Results



### Ordering Information

<b>Prod.#</b>	<b>Product description</b>
8479	Reveal 3-D for Total Milk Allergen

### Neogen Corporation

800/234-5333 • 517/372-9200 • E-mail: [foodsafety@neogen.com](mailto:foodsafety@neogen.com)

### Neogen Europe

+44 (0) 1292 525600 • E-mail: [info@neogeneurope.com](mailto:info@neogeneurope.com)